

Tek-MLS Installation & Registration

Rev. 042226

This document is provided to help users through the Tek-MLS installation and registration process. Tek-MLS is structured differently compared to previous versions of Tekran software. At first launch, the software requires a registration step. This step creates a software key unique to your PC and instrument. This process assures the software is registered properly for automatic notification of new software updates. These may be directly downloaded and installed (requires active internet connection).

To have access to the full functionality of the software, the software will need to have an active license. These may be purchased in one-year or multiple-year increments. If you are at the end of your license interval – the software will want the user with a popup notification that the license period has expired. Once the license expires, the software will have limited access to premium features.

PC Requirements:

Minimum System Requirements (For Single Tek-MLS Instance):


1. CPU: Intel i3 8th Generation or equivalent AMD CPU
2. OS: Windows 10 1809 or higher
3. Storage: Hard disk drive with enough storage spaces (100GB or higher)
4. Graphics Card: Integrated Graphics Card, GeForce 770, R9 280X equivalent or higher
5. Display: One monitor with 1366 X 768 resolution or higher
6. RAM: 8GB or higher

Recommended System Requirements (For up to 2 Tek-MLS Instances):

1. CPU: Intel i5 8th Generation or equivalent AMD CPU
2. OS: Windows 11 with latest updates
3. Storage: Solid-state drive with enough storage spaces (500GB or higher)
4. Graphics Card: Integrated Graphics Card, GeForce 770, R9 280X equivalent or higher
5. Display: One monitor with 1920 X 1080 resolution or higher
6. RAM: 16GB or higher


Installation Procedure:

1. Download Software – the Tek-MLS software may be downloaded from Tekran’s website at the hyper link listed below, or visiting the Tekran website downloads page.
2. Once downloaded, double-click on the file to install the Tek-MLS software on your PC.
3. Click the “Accept the Agreement” and hit “Next” button. Follow the on-screen prompts to install the software.

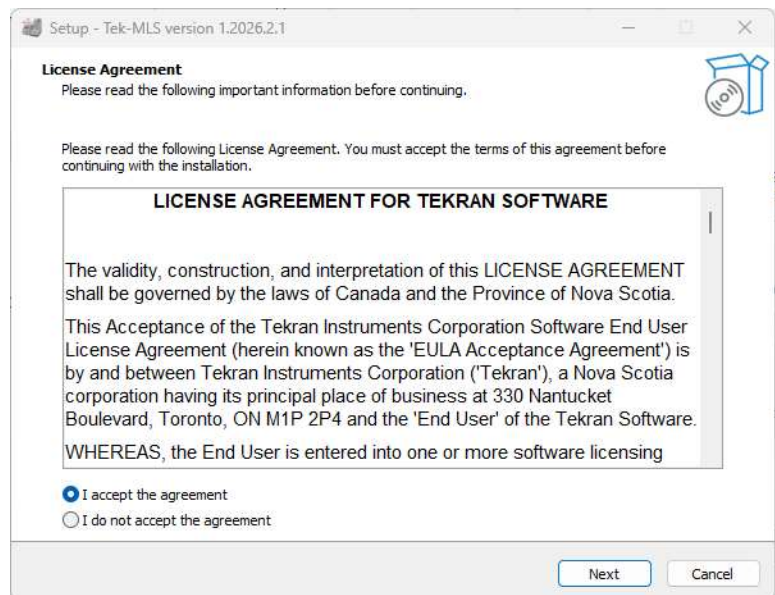

Software Download Link: https://www.tekran.com/files/Software/TekMLS_x64.exe

- Installing the software will require License Registration. Note this needs to be installed on the PC connected to the 2600 or 2700 instrument. The software will attempt to read the instrument serial number. If this is not possible, the registration will request manual entry of the instrument serial number. Select Instrument Type and enter Serial Number.

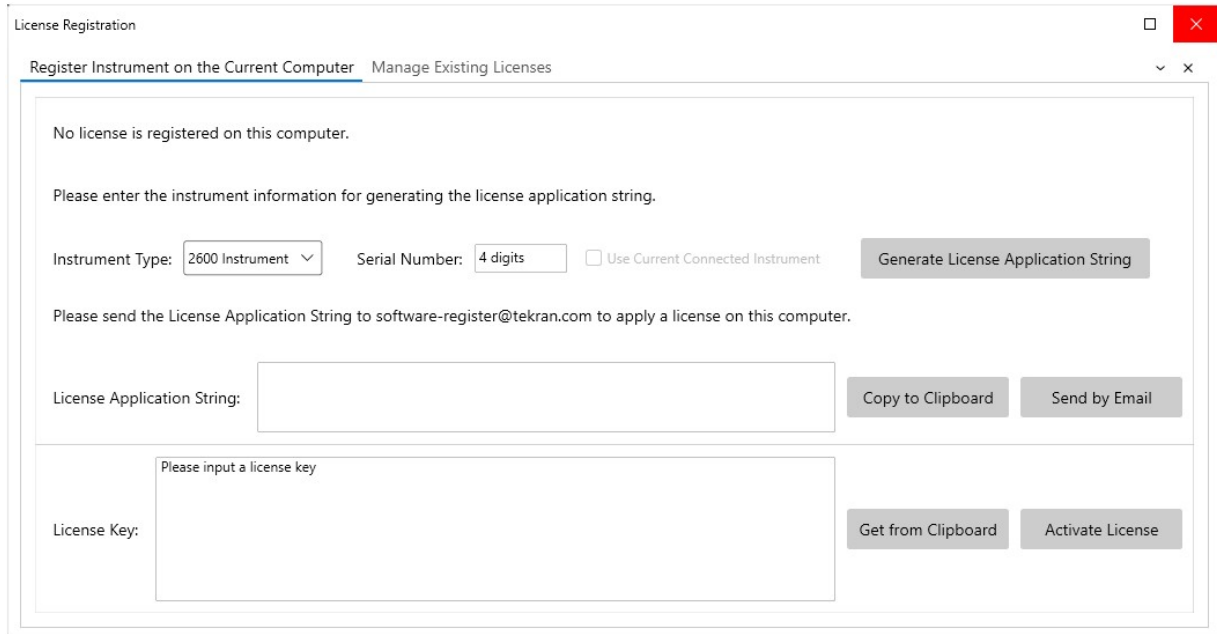

Warning: Tek-MLS is only compatible with the Tekran 2700 and 2600-IVS series instruments. Older 2600-STD systems with peristaltic pumps are not compatible with the Tek-MLS software.

- Once the instrument and serial number are entered, select the “Generate License Application String” button. This will create the application string in the box. Either hit the Send By email button or Copy to Clipboard.

Note: You may not use “Ctrl C” to copy the License Application String to the clipboard. The “Copy to Clipboard” button must be used for this function.

- After sending this License Application String to Tekran, it will be validated and a License Key will be sent back to you. Paste the License Key into the License Key box. Again, use the “Get from Clipboard” button for this function.
- Select the “Activate License” button. Congratulations on verifying your new Tek-MLS software! After the license period has ended, the software will provide an on-screen warning. There is a 30-day grace period to renew the license and maintain full software functionality. When trying to initiate a run with an expired license, a message similar to the below window will pop up.

